

MC-Media Pad® *E. coli* & Coliform test


Interpretation Guide

The MC-Media Pad® is a convenient method for indicator organism testing. Compliant with international food and beverages standards (AOAC-PTM, MicroVal, ISO 16140), MC-Media Pad® tests streamline your microbial quality control workflow.


Accurate results in 4 steps




1. Open the cover film diagonally

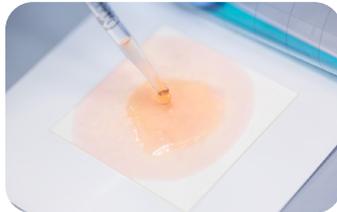

2. Inoculate your sample, it will diffuse automatically

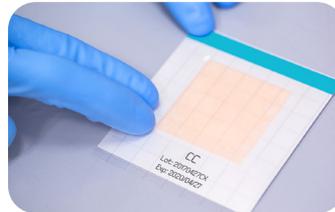

3. Close the cover film and incubate

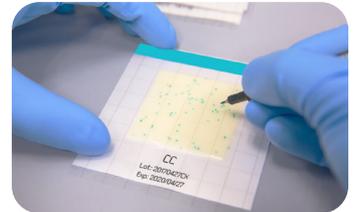


4. Count your results



Coliform

E. coli


Enterobacter aerogenes
ATCC 13048 and *Escherichia coli* ATCC 25922
50 CFU's

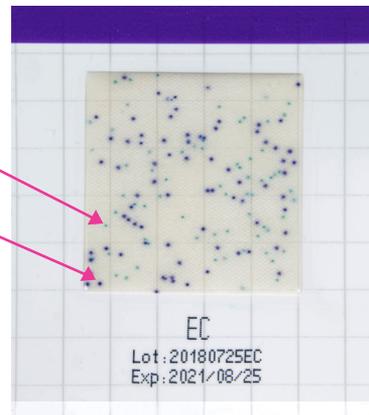

Enterobacter aerogenes
ATCC 13048 and *Escherichia coli* ATCC 25922
150 CFU's



Enterobacter aerogenes
ATCC 13048 and *Escherichia coli* ATCC 25922

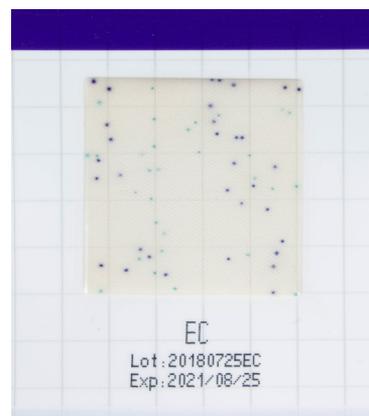

Enterobacter aerogenes
ATCC 13048 + *Escherichia coli* ATCC 25922 63 CFU's


With MC-Media Pad®, it is easy to see when a dilution is needed.



Dilution for easier read-out



Dilution factor 1:10

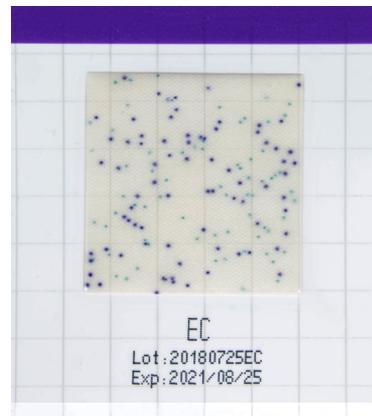

Dilution for easier read-out



Dilution factor 1:10



Ordering information	Product name	Packaging
1.32357.0002	MC-Media Pad® <i>E. coli</i> and Coliform	100 pcs

Please visit: SigmaAldrich.com/MC-MediaPad

